


NMR softwares installation



Mestrelab Mestrelab (EPFL site licences)

1. Download and install Mestrelab software for your operating system from Mestrelab [webpage](#)

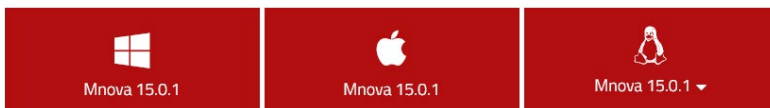

2. Download and unzip EPFL launcher from [this webpage](#) (need gaspar login on the EPFL network)

- [Launcher NMR PC Version](#)
- [Launcher NMR Mac Version](#)
- [Launcher NMR Linux Version](#)

3. Always start Mestrelab with the EPFL launcher (on the EPFL network or with VPN)




Topspin (Bruker) (Free academic licences)

1. Create a Bruker account with an academic address from [this webpage](#) (or sign in if you already have one)
2. Download Topspin latest version for your operating system from [this webpage](#) and install it (can take some time)
3. Request an academic license from [this webpage](#)

[REQUEST ACADEMIA LICENSE](#)

NMRium NMRium (Free)


1. Open [NMRium.org](#) [webpage](#)
2. Drag and drop your spectrum folder to open it

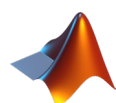

RBNMR + Matlab (Matlab licence needed)

1. Download RBNMR function (Nils Nyberg) from MatWorks [webpage](#)
2. Add rbnmr script to your Matlab functions folder
3. Run it with `X=rbnmr(datapath)`

NMR softwares installation

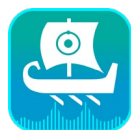

Jeol JASON (Licence need to be purchased)

1. Software and licence can be downloaded / bought on this [webpage](#)